

ReadiLink™ Protein Conjugation Stop Buffer

Catalog number: 5400
Unit size: 1 ml

Component	Storage	Amount
ReadiLink™ Protein Conjugation Stop Buffer	Refrigerated (2-8 °C)	1 vial (1 ml)

OVERVIEW

ReadiLink™ Protein Conjugation Stop Buffer is used to stop the protein conjugation reactions that use succinimidyl ester, sulfonyl chloride or isothiocyanate. It can be used directly without dilution.

AT A GLANCE

Important Warm up ReadiLink™ protein conjugation stop buffer at room temperature before use.

SAMPLE EXPERIMENTAL PROTOCOL

1. Prepare your protein (antibody) conjugation reaction as needed.
2. Stop Conjugation reaction by adding ReadiLink™ protein conjugation stop buffer at 1:50 ratio (such as for a 100 µL conjugation reaction, add 2 µL ReadiLink™ protein conjugation stop buffer).
3. Incubate at room temperature for 10 minutes.
4. The labeled protein (antibody) is now ready to use.

Note It is not necessary to purify conjugated proteins (antibodies) with a column after the conjugation reaction is stopped with ReadiLink™ protein conjugation stop buffer if the conjugation is used within 1 month.

Note For long term storage, it is recommended purifying the conjugated proteins (antibodies) with desalting column. And the protein conjugate should be stored at > 0.5 mg/mL in the presence of a carrier protein (e.g., 0.1% bovine serum albumin). For longer storage, the protein conjugates could be lyophilized or divided into single-used aliquots and stored at ≤ -20°C.

EXAMPLE DATA ANALYSIS AND FIGURES



Figure 1.
Mechanism for bioconjugation of dye to a peptide using succinimidyl ester.

DISCLAIMER

AAT Bioquest provides high-quality reagents and materials for research use only. For proper handling of potentially hazardous chemicals, please consult the Safety Data Sheet (SDS) provided for the product. Chemical analysis and/or reverse engineering of any kit or its components is strictly prohibited without written permission from AAT Bioquest. Please call 408-733-1055 or email info@aatbio.com if you have any questions.